

OMERO.web : First Steps

1 Login into OMERO.web

<https://omero.cellnanos.uni-osnabrueck.de/webclient/login/>




2 Navigate to data

Clicking on the group name in the top-left corner (next to your name) will reveal the groups your account belongs to



To see other people's data, select the name of the group you wish to browse and then the user inside that group whose data you wish to see

You can browse **folders** in the left-hand pane: Image folders are called **Datasets** and they are within **Projects**

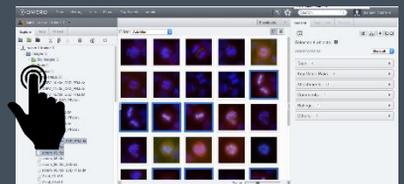

When a Dataset is selected, Image thumbnails are shown in the centre panel

3 How to open an image in OMER.iviewer

Double click on an image thumbnail or on the name will open the image in a viewer.

Viewing Images does not download the whole Image to the client. Only the viewed Image plane is rendered from the original Image file on the server and sent back to the OMERO.web client.

